

Most Affordable ATP Swab!

Ultrasnap™ ATP Swab

Order #US2020

Designed exclusively for the systemSURE II™ luminometer

Ultrasnap™ ATP swab is an integrated device using a unique liquid-stable reagent and Snap-Valve™ technology to deliver unbeatable accuracy and reproducibility. Simply swab an area, snap, squeeze and read the results with the systemSURE II™ luminometer in just 15 seconds. Ultrasnap™ swab and the systemSURE II™ puts the most accurate, reliable and affordable ATP-monitoring system in the palm of your hand.

Key Features

- Costs up to 50% less than competitor ATP swabs
- All-in-one reagent device
- Pre-wetted swab
- Tests both solid surfaces and liquid samples
- Unique liquid-stable reagent
- Unprecedented accuracy and reproducibility
- Robust, tolerant to temperature abuse
- 12-month shelf life at refrigerated temperature/
6-week shelf life at room temperature

User Benefits

- Significant cost savings
- Precise and sensitive
- Consistent and reliable
- Ready to use
- Simple, quick and convenient
- No reagent wastage
- Instant results within 15 seconds
- Objective measure of cleanliness
- Double your testing at no extra cost



941 Avenida Acaso, Camarillo, CA 93012
tel: 805.388.8007 ext. 300
e-mail: info@hygienausa.com
web: www.hygienausa.com

How to use:
solid surfaces or liquid samples

Snap




Squeeze



Insert device




Swab

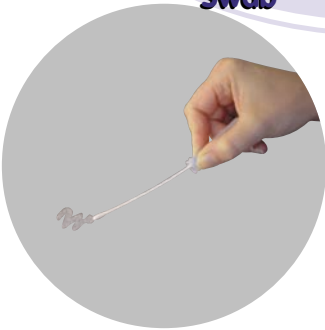

Use the Ultrasnap™ ATP swab and systemSURE™ luminometer to:

- Determine the cleaning efficiency and hygienic status of surfaces in food and beverage processing plants, restaurants, supermarkets and other industries
- Detect contamination in liquid samples particularly in water treatment plants or Clean In Place (CIP) systems
- Ensure food safety, improve product quality and reduce costs

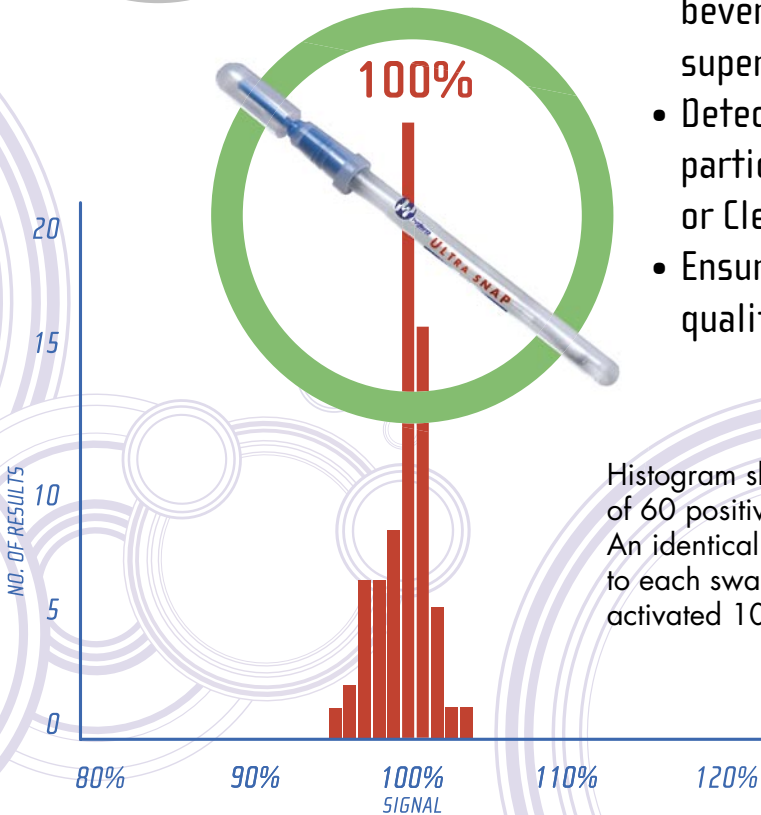

Histogram showing the reproducibility of 60 positive Ultrasnap™ ATP swab results. An identical amount of pure ATP was added to each swab tip; 6 different operators each activated 10 replicate devices.



hygiena

Rapid Solutions for Food Safety

941 Avenida Acaso, Camarillo, CA 93012

tel: 805.388.8007 ext. 300

e-mail: info@hygienausa.com

web: www.hygienausa.com